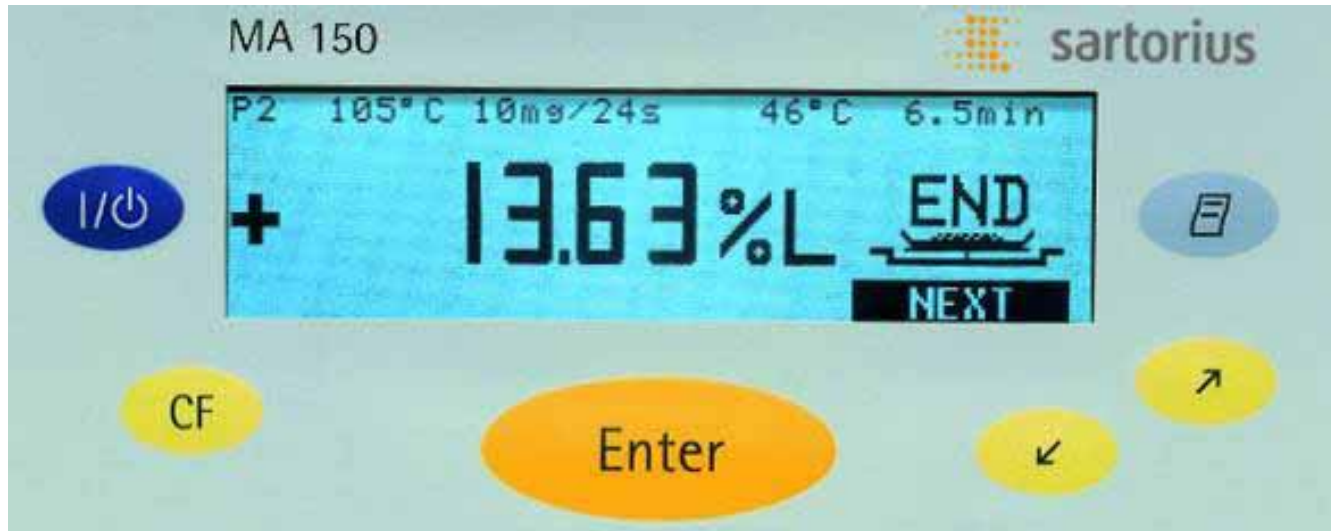
SARTORIUS MA150 的操作方法:

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| 步骤 | 按键 | 显示 |
|-----------------------|---|---|
| 1. 开机 |  |  |
| 2. 选择 PROGAM , |  | <u>PROGAM</u> |
| 3. 确认 PROGAM |  | P1 / P2 / P3 / P4 / P5 Heating Program Standby temperature Bar graph for weighing-in sample Start analysis End of analysis end Display mode Print intermediate results Factory setting |

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| | | |
|------------------------------|---|--------------------------|
| 1. 选择 Heating Program |  | O Heating Program |
| 2. 确认 Heating Program |  | O Standard drying |
| 3. 确认 Standard drying |  | Temperature 105°C |
| 4. 设定加热温度 (此例中为 130 度) | 重复按  | Temperature 1××°C |
| |   | Temperature 12×°C |

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| 步骤 | 按键 | 显示 |
|-------------|---|--|
| 8. 按 CF 键退出 |  | O Heating Program |
| 9. 按 CF 键退出 |  | Heating Program Standby temperature Bar graph for weighing-in sample Start analysis End of analysis end Display mode Print intermediate results Factory setting |

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O Bar graph for weighing-in sample

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Heating Program
Standby temperature
Bar graph for weighing-in sample
Start analysis
End of analysis end
Display mode
Print intermediate results
Factory setting
显示

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Heating Program
Standby temperature
Bar graph for weighing-in sample
Start analysis
End of analysis end
Display mode
Print intermediate results
Factory setting

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O Start analysis

2. 确认 **Start analysis**

With stability ,after Enter ket
W/O stability , after closing

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| | | |
|-------------|---|--|
| 9. 按 CF 键退出 |  | Heating Program Standby temperature Bar graph for weighing-in sample Start analysis End of analysis end Display mode Print intermediate results Factory setting |
|-------------|---|--|

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